

EYE WASH BOTTLE

An eye wash bottle is a portable device filled with sterile saline solution, used to quickly flush out irritants, chemicals, or debris from the eyes in emergencies.



ABOUT US

Waltzer India is a trusted name in industrial safety, offering high-quality eye wash stations, safety showers, PPE, and custom solutions. Serving industries like pharma, chemical, and manufacturing, we combine innovation with global safety standards to protect people and workplaces.



Scan here

CONTACT

Waltzer India



A-Balaji Market, Hawa Bangala
Indore- 452009 Madhya Pradesh



Waltzify.com



+91 95225 82422
sales@waltzerindia.com

EYE WASH SOLUTION & BOTTLE



Simple. Safe. Effective

EYE WASH SOLUTION

Formulated for immediate eye irrigation in emergency situations involving dust, chemicals, or foreign particles. This 500 ml sterile solution is designed to flush contaminants and soothe the eye area without causing irritation.



How to Use

1. Fill the bottle with sterile eye wash solution.
2. Attach the nozzle and make sure it's secure.
3. Tilt your head toward the affected eye over a sink.
4. Place nozzle near the eye and squeeze the bottle gently.
5. Keep eye open and blink to flush out irritants.
6. Flush for 10-15 mins if exposed to chemicals.

Important Notes:

- For external use only
- Do not reuse – single-use product
- Keep away from children
- Store in a cool, dry place



Why Use Eye Wash Solution?

Because accidents don't wait – and your eyes don't get a second chance.

- Immediate relief
- First-Aid essential
- Sterile & Safe
- Prevents permanent damage

